

Lentiviral Transduction of CRC Organoids and CAFs

Contact address: Henner F. Farin (Georg-Speyer-Haus, Institute for Tumor Biology and Experimental Therapy, Frankfurt/Germany; email: h.farin@georg-speyer-haus.de)

Version 1.0 Version Date: 11 April 2023, generated by Benardina Ndreshkjana

Introduction

Lentiviral transduction is an efficient method to introduce transgenes into mammalian cells for robust and stable expression. The possibility to express reporter proteins in organoids and cancer-associated fibroblasts (CAFs) using lentiviral vectors has increased the opportunities for the design of patient-derived cell assays. Handling of lentiviral constructs requires appropriate permissions and laboratory infrastructures to work under *Biosafety level 2* conditions.

Purpose

This standard operating procedure (SOP) describes the procedures for lentiviral transduction of CRC organoids and CAFs. Virus production and titration is performed as described in *Michels et al., 2020, Cell Stem Cell 26, 1–11*.

1. Equipment, materials and chemicals

1.1 Equipment

- Cell culture incubator adjusted at 37 °C and 5 % CO₂
- Centrifuge (Rotanta 460R, Hettich)
- Evos Fl microscope with 2X, 4X, 10X objectives
- Pipettes with sterile tips
- Pump and aspiration pipettes
- Water bath or heating block
- Pipette boy

1.2 Materials

- 50-ml Falcons
- 5 ml Eppendorf tubes
- Sterile 6 cm and 10 cm cell culture dishes
- Sterile 12-well cell culture suspension plates (Greiner, #665102)
- 5- and 10-ml pipets
- Parafilm M

1.3 Chemicals

- Matrigel™ (Corning, #356231)
- Y-27632 Rho kinase inhibitor (Biotrend, #Y5301)
- PBS (Thermo Fisher, # 10010015)
- 0.1 % gelatin in PBS (Sigma-Aldrich, #G2500-100G)
- Trypsin (Thermo Fisher, #12605-010)
- Polybrene Transfection Reagent (Sigma-Aldrich, #TR-1003-G)
- Puromycin (InvivoGen, #ant-pr-1)
- Accutase (Life Technologies, #A1110501)

2. Medium recipes

2.1 Washing medium (WM)

Component	Concentration (stock)	Amount
DMEM (Thermo Fisher, #31966-047)	1x	500 ml
Bovine Albumin Fraction (Thermo Fisher, #15260-037)	7.5%	2.65 ml
Pent/Strep (Thermo Fisher, #15140-122)	100x	5 ml

2.2 Advanced DMEM+++

Component	Concentration (stock)	Amount
Advanced DMEM/F12 (Thermo Fisher, #12634-028)	1x	500 ml
Glutamax (Thermo Fisher, #35050038)	100x	5 ml
HEPES (Thermo Fisher, #15630056)	1M (100x)	5 ml
Pent/Strep (Thermo Fisher, #15140-122)	100x	5 ml

2.3 Medium for tumor organoid expansion

Component	Concentration (stock)	Amount
Advanced DMEM +++	1x	35 ml
*R-spondin (in house prepared CM)	1x	10 ml, 20% (v/v)
**Noggin (in house prepared CM)	1x	5 ml, 10% (v/v)
B-27 (Thermo Fisher, # 12587010)	50x	1 ml
N-acetylcysteine (Sigma-Aldrich, #A9165)	5M	125 μ l
Epidermal growth factor (Peprotech, #AF-100-15)	500 μ g/ml	5 μ l
ALK5 inhibitor A83-01 (Tocris, #2939/10)	500 μ M	50 μ l
p38 inhibitor SB202190 (Sigma-Aldrich, #S7076)	30 mM	16.6 μ l

*R-spondin1-conditioned medium was obtained from R-spondin producing cells (293T-HA-Rspol-Fc, from Calvin Kuo lab) and used under approved MTA.

**Noggin conditioned medium was obtained from Noggin producing cells (HEK293-mNoggin-Fc, from Hans Clevers lab) and used under approved MTA.

2.3 Medium for tumor fibroblast expansion

Component	Concentration (stock)	Amount
DMEM (Thermo Fisher, #31966-047)	1x	500 ml
Fetal Bovine Serum (Sigma-Aldrich, #F7524)	10x	60 ml
Pent/Strep (Thermo Fisher, #15140-122)	100x	5ml

3. Transduction of organoids

- Thaw the virus (200 μ l/transduction) and Matrigel on ice. Use 1 full well of organoids that was passaged at least 5 days before transduction.
- Remove the medium, add 1 ml of cold washing medium (WM) and resuspend the Matrigel drops using a 1000 μ l pipette.
- Collect the organoids in a 5 ml Eppendorf tube.
- Mechanically shear the organoids using a 10 μ l tip on top of a 1000 μ l tip.
- Vigorously pipette up and down 20-30 times.
- Fill up with cold WM and spin down at 1200 rpm for 5 min at 4 °C.
- Remove the supernatant of the pelleted organoids.
- Add 1 mL of Accutase and pipette up and down a few times.
- Incubate for 3 min at 37 °C in a water bath or heating block.
- Pipette up and down and check under the microscope if most cells are singularized.

- Fill up the tube with WM and pellet the organoids at 1500 rpm for 5 min at 4°C.
- Remove the supernatant and resuspend the cell pellet in 1 ml of Advanced DMEM+++ medium (adv DMEM).
- Divide the organoid suspension in 750 µl for the transduction and 250 µl for the control in a 12 well suspension plate.
- To the transduction add 200 µl of virus and 50 µl of adv DMEM medium.
- To the control add 750 µl of adv DMEM medium.
- In both wells add 1:1000 Y-27632 Rho kinase inhibitor and 1:1000 Polybrene.
- Wrap the plate with parafilm and centrifuge at 2200 rpm for 1 h at 32 °C.
- Remove the parafilm and place the plate in the incubator for 3 h.
- Collect each well in a 5 ml Eppendorf tube with 5 ml of WM and centrifuge at 1500 rpm for 5 min at 4 °C.
- Repeat the washing step 2 times.
- Remove the supernatant and add 20 µl of adv DMEM medium.
- Resuspend the transduction mix with 80 µl Matrigel and make 8 drops of 10 µl (4 drops/well) in a 12 well suspension plate.
- Resuspend the control in 40 µl Matrigel and make 4 drops of 10 µl (2 drops/well). In one well add Puromycin and in the other not.
- Place the plate in the incubator for 20-30 min.
- Add 1 ml of pre-warmed organoid medium, supplemented with 1:1000 Y-27632 Rho kinase inhibitor/well and place it back in the incubator.
- After 3 days start selection with 1:1000 Puromycin (leave one control well without Puromycin).
- Usually, the organoids are ready for passaging after 1-2 weeks.

4. Transduction of fibroblasts

- Thaw the virus on ice (200 µl/transduction). Usually use one full 10 cm culture dish with fibroblasts for 5 transduction reactions.
- Add 1 ml of 0.1 % gelatin in PBS to five 6 cm culture dishes and let them solidify for 20-30 min in the incubator.
- Remove the medium from the fibroblasts, add 3 ml of PBS and remove it as well.

- Add 1 ml of Trypsin and place it back in the incubator for 4-5 min.
- After 4 min check under the microscope if the cells started to detach and collect with 4 ml of fibroblast medium in a 5 ml Eppendorf tube.
- Centrifuge at 1500 rpm for 5 min at 4 °C and remove the supernatant.
- Collect the cell pellet in 1 ml fibroblast medium.
- Take up the gelatin from the 6 cm dishes and add 3.6 ml of pre-warmed fibroblast medium with 1:1000 Polybrene.
- In each plate add 200µl of the fibroblasts and 200 µl of virus and place them back in the incubator.
- After 5-6 h remove the medium and add 4 ml of fibroblast medium (check before if fibroblasts are attached).
- After 3 days start selection with 1:4000 Puromycin.
- When the plates are confluent, split the 5 plates in 2 of 10 cm gelatin-coated dishes. After each passage slowly increase the Puromycin concentration to 1:2000.